Central Research Laboratory

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Guidance for New Users FY2026
Research Equipment Section
Central Research Laboratory

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1) Overview: Usage of Research Equipment Section,
CRL

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User registration process at Research Equipment Section, CRL

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For on site attendee

1. Attend this lecture
2. Download "Application Form" and submit it by e-mail attachment (hqcr@belle.shiga-med.ac.jp)

For Zoom on line attendee

1. Attend this online Zoom guideline lecture
2. We will provide a unique link in the Zoom chat during the lecture
3. Go to the link and download both "Survey form" and "Application Form"
4. Fill out both forms and send them by e-mail attachment for approval

* By annual renew application, the user registration is maintained.

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Why is the user registration required?

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1. Request from National University Corporations

- ・Compliance with [Industrial Safety and Health Act](#)
- ・[Experimental safety management and education](#)

2. Stricter regulations in life science research

- ・Recombinant DNA experiments: Act on the Conservation and Sustainable Use of Biological Diversity through Regulations on the Use of Living Modified Organisms, etc.
- ・Experimental animals: Act on Welfare and Management of Animals, etc.
- ・Experiments handling pathogenic microorganisms: Act on the Prevention of Infectious Diseases and Medical Care for Patients with Infectious Diseases, etc.

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Request in the use of CRL facilities

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<Basic Rules>

1. Fill out the booking sheet of each instrument that you plan to use; Write the proposed time of use, your name and belonging, and the extension number.
2. Prepare any consumables, tools, chemicals, etc. for your own use.
3. Strictly follow the operating rules to use instruments and analyzers.
4. Fill out the user record sheet after every use (This is important for calculating the user charge).
5. Maintain the facility in a clean condition after use.

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Cautionary notice for booking equipment

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1. Write the extension number on the booking sheet.

Users should communicate with other users and the CRL staffs.

2. Do not make booking for inappropriately long period of time.

The equipment booking may be canceled if the equipment is not used after 30 minutes of proposed starting time, especially for equipment that is commonly used by many users.

3. Follow the booking rules for each equipment.

4. Follow the CRL rules and strictly observe the start/end time for use.

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Follow the CRL operating rules to use equipment

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- ・When using equipment at CRL for the first time, **any users MUST ask the CRL staffs** how to operate it.
- ・Online operation manuals are available on the CRL website (http://www.crl.shiga-med.ac.jp/home/kiki_bumon/g_book/contents.html). (Sorry, we will create the English version soon)

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1. The CRL facility is open 24 hours a day.
For overtime and holiday use of the locked laboratories, contact the CRL staff in advance to get the key for the labs.
2. There are laboratories statutorily designated as controlled areas: Genetic Engineering Lab, Pathogenic Virus Lab, Animal X-ray Fluoroscopy Lab.
・ Permission and registration are necessary to enter the above labs.
・ **Strictly follow the regulations defined by SUMS and CRL.**
3. CRL provides "Special service" for several instruments and analyzers.
Specific equipment is operated only by the CRL staffs for the support of researchers. To analyze the samples by using the equipment, bring the samples to the staff with completed application form.

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- ・Central Research Laboratory Seminar (TBA)
- ・Technical Seminar (TBA)
- ・Intensive Course Seminar (Sep 8-11, 2026)
in "Basic Science Fundamentals & Multidisciplinary Seminars"
Participants in the above seminars can obtain credits for the PhD course lectures.
Details are indicated at the CRL website.
(<http://www.crl.shiga-med.ac.jp/home/seminar/seminar.html>)

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Tips for users of Research Equipment Section, CRL

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1. Please visit the CRL website for better usage of Research Equipment Section
(<http://www.crl.shiga-med.ac.jp/index.html>)
2. Good communication with CRL staffs is very helpful for both users and staffs.

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On site attendee

After this lecture, download the application form and submit it to CRL via e-mail

On line Zoom attendee

We will display a link for download site of two forms to be submitted ("survey form" and "application form for CRL User Registration") in the Zoom chat box during this course. By sending them using email, your registration will be completed.

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SDS provide procedures for handling or working with that substance in a safe manner.

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Formaldehyde should be used under the fume hood in the CRL.
 Users should sign in a Formaldehyde Log after use.

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Evacuate from the fire zone if you think it dangerous.

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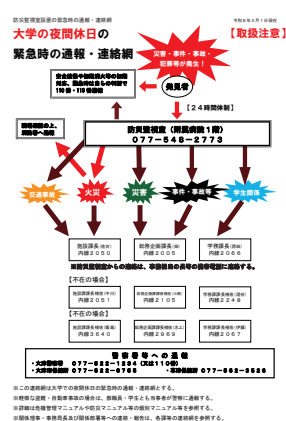
How to use the "recombinant DNA room" and
"pathogenic virus room"

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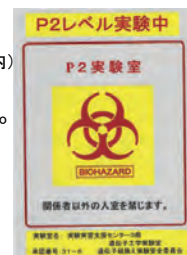
Act on the Conservation and Sustainable Use of Biological Diversity through Regulations on the Use of Living Modified Organisms(LMOs), so-called **Cartagena Law**

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Containment measures (Type 2 Use of LMOs)

Possible uses are those with containment measures to prevent to dispersal of LMOs in the environment (Uses in laboratories etc).

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Disposed E. coli solution without autoclaving.
Serious violation of Cartagena Law.

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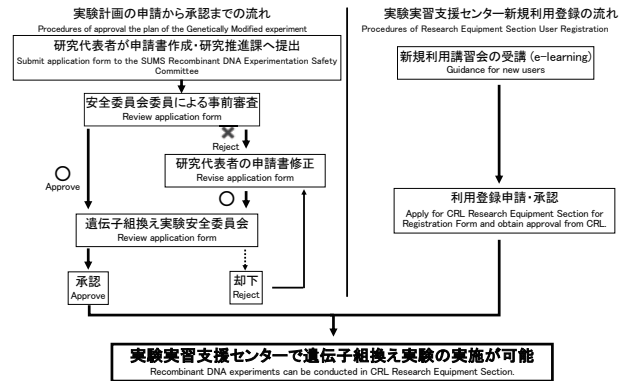
The government may order suspension of research.
Offender shall be punished by imprisonment or a fine, or a combination of these two.

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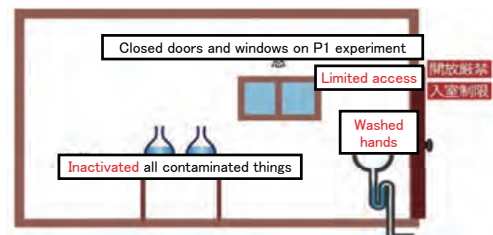
P1レベル拡散防止措置 (抜粋) Containment Measures of P1 Level Laboratory

P1Aレベル拡散防止措置 (抜粋) Containment Measures of P1A(experiment with Animal) Level Laboratory

遺伝子組換え生物等が逃亡その他拡散しない構造の容器に入れること
When a LMO is taken out of the laboratory in process of an experiment, the LMO shall be put in a container of the structure that prevents it from leaking or other dispersion.

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When using the equipment for the first time, please ask the CRL staffs how to use it.

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Do not hesitate to ask any questions to CRL staff.

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Do not leave the waste after autoclaving.

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

ウイルス実験室3
Pathogenic Virus Lab 3
(for Retrovirus, Lentivirus)

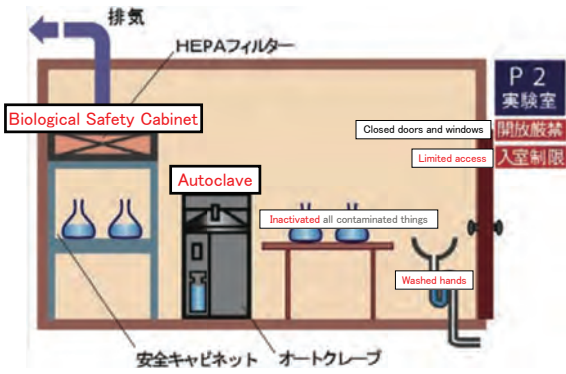P2レベル拡散防止措置(抜粋)
Containment Measures of P2 Level Laboratory

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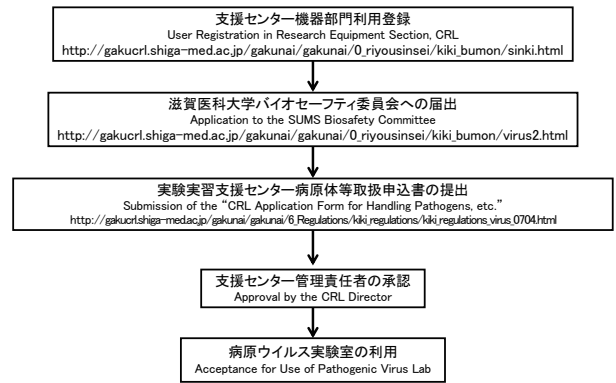4) How to start animal experiments at CRL

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Need to complete following:
1. Completed the training program in animal experiment.
2. Pass the qualification exam of animal experiment.
3. Obtain the certification on animal experiment.

動物生命科学センター登録申請
Apply for the registration to Research Center for Animal Life Science.

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Animal experiments in the CRL rooms

「動物の愛護及び管理に関する法律」、「滋賀医科大学動物実験規程」に従う。
Perform animal experiments under the Law Concerning the Protection and Control of Animals and the regulations of Research Center for Animal Life Science in SUMS.

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Approved rooms for animal experiments at CRL

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Technical staff, Central Research Laboratory

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You have 2 choices

- 1) prepare sequencing reaction and submit the sample for running
- 2) submit a plasmid for DNA sequencing.

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Download and fill in the Excel order form

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Submit samples with the order form to Room 307

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CRL's staff will run your samples and return the data

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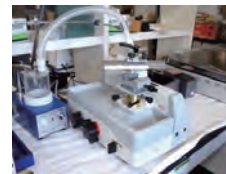
Yamamoto (Ext. 2304)
4F Room405

その他サービス
Additional support provided by CRL

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10x Genomics Single cell RNA-sequencing reagentの代理購入
Provide 10x Genomics Single cell RNA-sequencing reagent

qPCR plateの販売
Provide qPCR plates

MOLSIS Molecular Simulation & Informatics Systems Software

FlowJo software for Flow cytometer and FACS analysis

Leica LAS-X software for image processing

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A young researcher in the SUMS hospital wanted to determine gene mutations in the patient sample. The CRL staff helped designing primers, PCR, and DNA sequencing. The mutation was successfully detected.

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We are teaching how to do Western blotting.

Please ask us if you would like to know how to perform experiments. We may not help your issues directly, but will help your research as much as we can.

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X-ray Irradiator AcroBio RS320
1st floor of Research Center for Animal Life Science

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X-ray Diffractometer
1st floor of Central Research Laboratory

No risk of x-ray leaking from these generators

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X線照射装置使用の登録完了
Registration is completed for users of X-ray irradiators
(No need blood test and wearing a film badge)

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C-arm X-ray Fluoroscope ARCADIS Avantic Gen2
1st floor of Central Research Laboratory

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学内登録申請(まとめ)
Registration as a radiation worker in SUMS

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担当: 山内(内線2304)
Technical staff: T. Yamamoto (Ext. 2304),

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Introduction of each equipment available at CRL

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
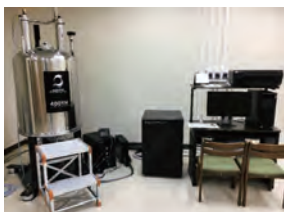
NMR can determine the molecular structure of a sample placed in a strong magnetic field by nuclear magnetic resonance.

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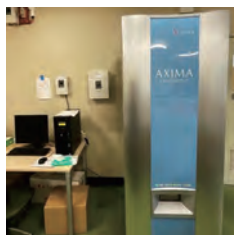The device is capable of measuring molecular weight of small molecules.

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This device can measure molecular weight of gas molecules.

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HPLC can separate and quantify components dissolved in a liquid sample.

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Compatible to 6-384 well Plates for measuring absorption, fluorescence, and luminescence.

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Nucleic acid, protein, etc. can be quantified with a small sample of 1 - 2 μ L.

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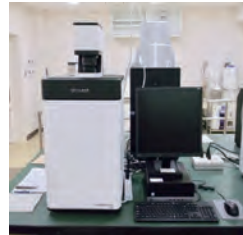
Ask the CRL staff which rotor will be compatible.

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
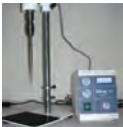
CRL 3rd floor

Various Cell Homogenization Equipment

Ultrasonic Cell Crusher

Homogenization with beads Minilys

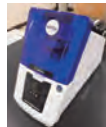Polytron Homogenizer

Vacuum Freeze Dryer

CRL 4th floor

Confocal Laser Scanning Microscope

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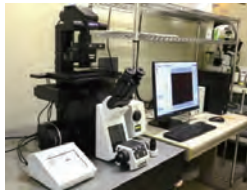
Confocal laser scanning microscopy is a technique for obtaining high-resolution optical images with depth selectivity.
Please ask CRL staff which microscope will be suitable for your experiment.

CRL 4th floor

Other microscopes

1. Inverted Fluorescence Microscope IX83 (Olympus)
2. DeltaVision High-resolution Fluorescence Microscope (GE)
3. All-in-one Fluorescence Microscope BZ9000 (Keyence)
4. Eclipse Ni-E (Nikon)

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
Ask us which microscope will be suitable for your experiment.

CRL 4th floor

New

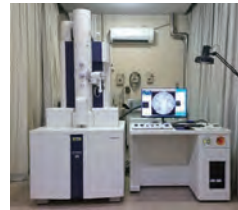Transmission Electron Microscope HT7800 (Hitachi)

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CRL provides sample preparation service for observation of the microstructure of biological specimens with an electron microscope.

Scanning Electron Microscope JSM-7505FA (JEOL) JSM-6010LA (JEOL)

CRL 4th floor

Ultramicrotome ULTRACUT E (Reichert-Jung)

Making thin sections for electron microscope

Draft Chamber

Use organic solvent under the draft chamber

CRL 4th floor

Automated Tissue Processor
Paraffin embedding

Tissue Embedding Console
Making paraffin blocks

Microtome
Making paraffin sections

Cryostat
Making frozen sections

CRL 4th floor

New

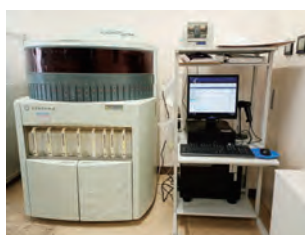
Digital Slide Scanner NanoZoomer S60v2

Virtual slide scanner

New

Automated Immunohistochemical Staining System VENTANA DISCOVERY ULTRA

Automate immunostaining

CRL 4th floor

Flow Cytometer

1. LSRFortessa X-20 (BD)
2. FACSCanto II (BD)
3. CytoFLEX S (Beckman)
4. FACS Calibur (BD)
5. Attune (AB)

Cell analysis

FACS

1. FACS Aria Fusion 5-Laser (BD)
2. FACS Aria Fusion 4-Laser (BD)

Cell analysis and sorting

Ask us which equipment will be suitable for your experiment.

CRL 4th floor

**Magnetic Cell Sorter
autoMACS (Miltenyi Biotec)**

Cell separation using antibodies

**Suspension Array System
Bio-Plex200 (Bio-Rad)**

Measuring cytokines using special beads

**Extracellular Flux Analyzer XFe24
(Seahorse Bioscience)**

Measuring metabolism in cultured cells

CRL 4th floor

CO₂ Incubator and Clean Bench

Users may request use of CO₂ incubator or clean bench for cell culture.

CRL 4th floor

**Large Capacity Liquid Nitrogen Cryopreservation
Equipment (MVE Heco1536P)**

Store cells in liquid nitrogen gas

Liquid Nitrogen

Users can buy liquid nitrogen from CRL

Milli-Q and Elix Water System (2nd and 4th floor)

Users can buy pure water from CRL

RCALS 3rd floor

**in vivo Imaging System
NEWTON 7.0**

Detection of bioluminescence and fluorescence in live animals

**High-resolution Ultrasonic Echo
Imaging System**

Visualize *in vivo* anatomical, functional, and physiological data in real time using ultrasound echo.

**Tetrapolar Bioimpedance Spectroscopy
Analyzer ImpediVET**

Measurement of fluid status and tissue composition in live animals

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8) How to avoid unnecessary accidents?

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The rotor came off from the spindle and rolled around in the chamber. The rotor and spindle were severely damaged.

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The bucket came off from the running rotor at 40,000 rpm and crashed onto the chamber wall at a velocity of over 1,600 km/h. Relative Centrifugal Field is 193,000 g.

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The lid of the refrigerated centrifuge was left closed for an extended period after use. The lid became sticky and stuck shut due to frost and organic solvents volatilized from the samples.

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Leica TCS SP8 X
Confocal Laser Scanning Microscope

In 2023, someone broke the objective lens.
It cost 700,000 yen for replacement.

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Nikon/AX Confocal Microscope

The user wanted to see the edge of the slide and moved the objective lens upward to adjust focus. Then, the objective lens hit and bent the stage holder.

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-
1. If you cause any accident, immediately inform it to the CRL staff.
 2. If you notice anything unusual such as a trouble with equipment, immediately inform it to the CRL staff.
 3. If you are not sure how to use the equipment, ask the CRL staff about it. DO NOT use the equipment only on your own idea and judgment.
 4. In using the CRL facilities, the rules at CRL are prior to those at your lab.
 5. DO NOT follow the user who acts in incorrect manner and method.

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